tecnosoft SOFTWARE INSTALLATION GUIDE

Step 1 - Download TecnoUtility

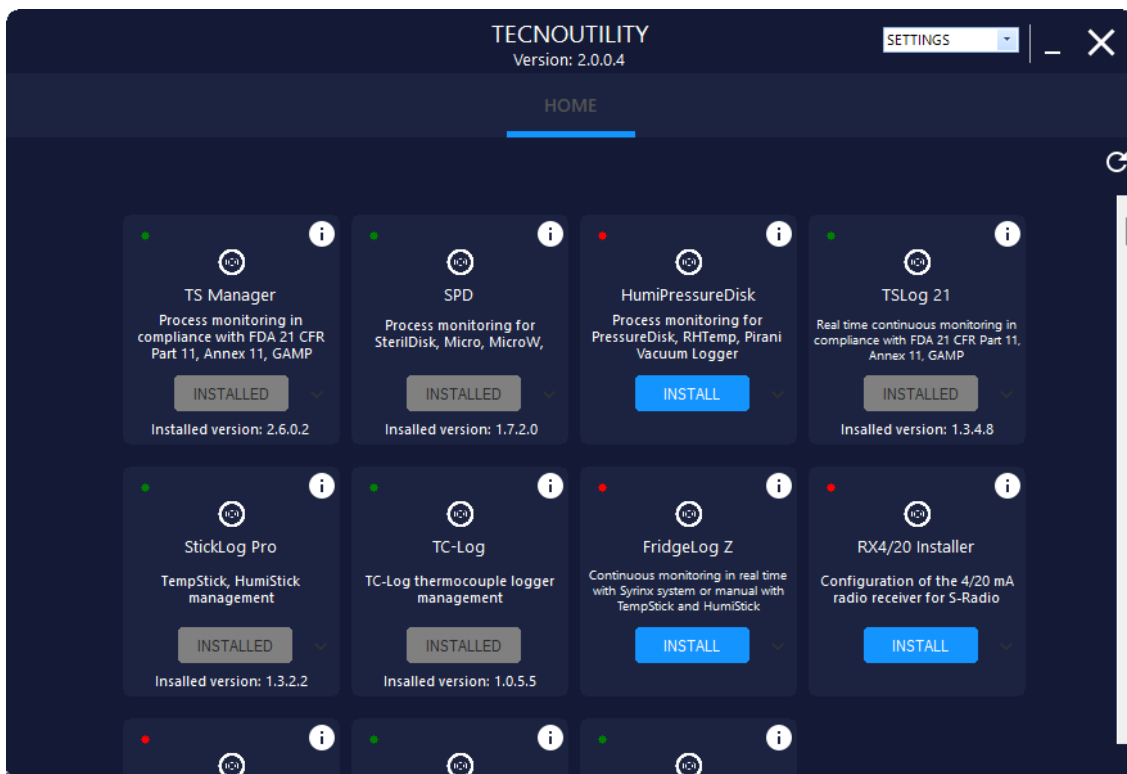
EN: <https://www.tecnosoft.eu/en/products/software-app/>
IT: <https://www.tecnosoft.eu/it/prodotti/software-app/>

Step 2 - Launch TecnoUtility

TecnoUtility is the all in one Tecnosoft app to manage all your monitoring software. Save it on your desktop and use it to access manuals, general documentation, video and software installation. When a new version is made available it will be prompted and you can upgrade it with a click.

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Step 3 - Install your software

Click on **INSTALL** under the desired software. If software is already installed and a new version is available, you can **UPDATE** it. Click on **i** for additional information.

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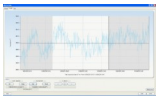
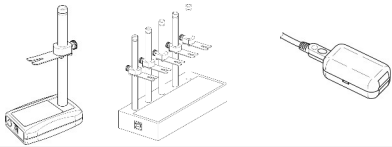
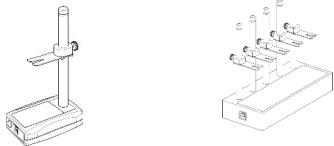
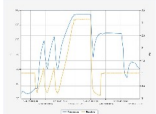
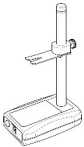
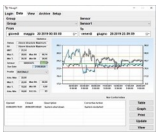
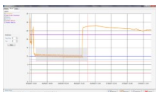
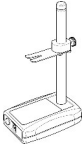
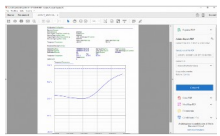
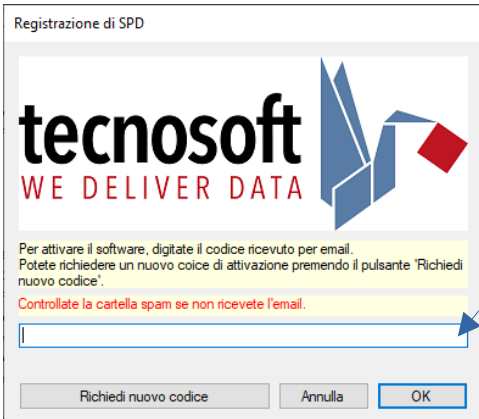
Software Table / Lista dei Software

Check the software you need to install according to the devices you have. Software with the  symbol require a USB protection key to be started and a full licence purchased.

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Step 4 - Registration

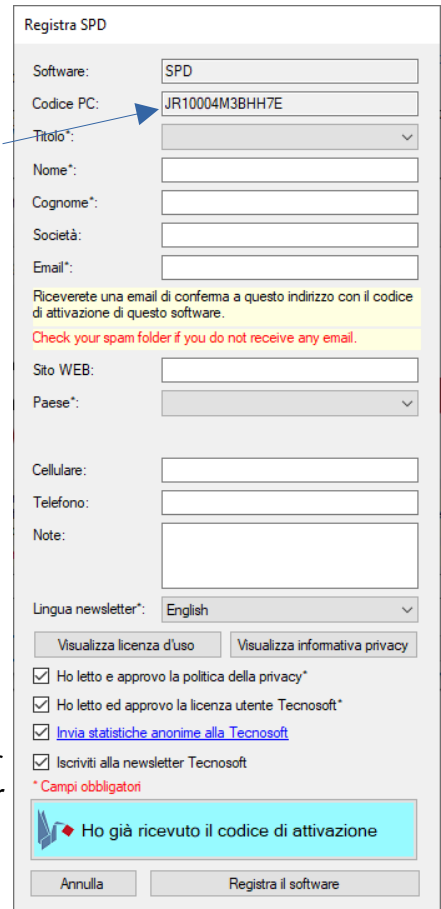
Registration is very simple and quick. Fill the form, send it, receive the activation code via email and start using your software!

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Your data here
I tuoi dati qui

Why should I have to register?

Registering the software is necessary in order to use the software: an activation code is sent by email with which you can activate your license. All collected data are stored following the European privacy directive (GDPR). Registration allows Tecnosoft to provide a more precise and immediate assistance service and to improve the software and prevent possible problems. Tecnosoft thanks you for your precious collaboration.

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CALIBRATION CERTIFICATES

Calibration certificates are available online only on our Support website.

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Useful Links

Support website: <https://support.tecnosoft.eu/en>
Registration FAQ:
<https://www.tecnosoft.eu/en/faqs/tecnosoft-support-website-registration/>
Certificates FAQ:
<https://www.tecnosoft.eu/en/faqs/#faq-category-certificates>

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For more Info and Support

Check our FAQs page:
<https://www.tecnosoft.eu/en/faqs/>
Especially:
<https://www.tecnosoft.eu/en/faqs/technical-assistance-service/>
or open a ticket on our Support website:
<https://support.tecnosoft.eu/en/assistance>

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TECNOSOFT SUPPORT WEBSITE

Access support.tecnosoft.eu to open a full range of support and assistance possibilities.

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You can download the reference certificate as well.

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Order the **PLATINUM** assistance for full coverage.

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Device

Problem

Lookup your model using the serial number

Can't find your product? Browse the catalog.

Question

UPLOAD ATTACHMENT

The file must not be over 2MB in size. Only jpg, png, gif, pdf and doc files are allowed.

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1 Select the intervention needed or the problem found

Request

- Standard calibration
- Extra calibration points
- Battery replacement
- Check

Problem

- Does not connect
- Wrong readings
- Mechanically broken
- Not calibrated
- No data recorded
- E1
- E2
- E3
- E4
- E5
- E6

2 List the serial numbers and add other notes

Find serial

Serial #

Device

Note

I don't know the serial number or the device name